

Green-Z® Spill Control Solidifier

(Chlorine-Free Formula)



At times, workers may be unintentionally exposed to blood and bodily fluid spills along with other harmful substances such as toxic chemicals, biological hazards, and chemotherapy agents. To help reduce those exposure risks, rely on Green-Z® for improved safety in your spill clean-up operations.

- Contains fast-acting, superabsorbent polymers
- Chlorine-free (safe for use on chemotherapy, glutaraldehyde and mercury spills)
- Eliminates splashing by solidifying and turning spills into semi-solid (gel) mass
- Cost-effective clean-up solution
- Single-Use Pouches Bottles

| SPECIFICATIONS | |
|----------------|---|
| Order No. | Description Case |
| 42019 | Green-Z® 21g single-use pouch 100/case |
| 42010 | Green-Z® 5 oz. shaker top bottle 24/case |
| 42012 | Green-Z® 15 oz. shaker top bottle 12/case |



Alky Solidifier

Rapidly Neutralizes & solidifies high pH spills

- Ammonium Hydroxide
- Potassium Hydroxide
- Sodium Hydroxide



Sprinkle Alky directly onto the spill, covering it completely. The spill will turn blue, then yellow when it is safely neutralized. Once it becomes a semi-solid gel, it is ready for clean-up with a scoop and scraper.

| SPECIFICATIONS | |
|----------------|--|
| Item No. | Description |
| BNP142 | ALKALI NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 142G |
| BNP425 | ALKALI NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 425G |
| BNP425/12 | ALKALI NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 425G BOX OF 12 |

Acid Lock Solidifier

Rapidly neutralizes & solidifies low pH spills

- Acetic Acid
- Nitric Acid
- Hydrochloric Acid
- Sulfuric Acid



Sprinkle Acid Lock directly onto the spill, covering it completely. The spill will turn pink, then yellow when it is safely neutralised. Once it becomes a semi-solid gel, it is ready for a cleanup with a scoop & scraper.

| SPECIFICATIONS | |
|----------------|--|
| Item No. | Description |
| ANP142 | ACID NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 142G |
| ANP425 | ACID NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 425G |
| ANP425/12 | ACID NEUTRALISING ABSORBENT POLYMER SHAKER TOP BOTTLE 425G BOX OF 12 |

